

Chemical resistance guide for Finite HD

It is important to note that the performance of these gloves may be different if mixtures of chemicals are involved.

Chemical	EN374 performance level			
Acetic acid 50%	2			
Acetone	0			
Acetonitrile	0			
Carbon disulphide	0			
Dichloromethane	0			
Diethylamine	0			
Heptane	6			
Hydrochloric acid 32%	6			
Methanol	0			
Nitric acid 40%	4			
Phosphoric acid 85%	6			
Sodium hydroxide 40%	6			
Sodium hypochlorite 13%	6			
Sulphuric acid 40%	6			
Sulphuric acid 98%	1			
Toluene	0			
Xylene	0			

EN374 Performance Levels

Performance Level	1	2	3	4	5	6
Breakthrough time (mins)	>10	>30	>60	>120	>240	>480

The information contained here is advisory and is intended to guide and inform only. This information is based on the view of BM Polyco Ltd in light of all available data and is not subject to any guarantee on our part regarding the performance of a particular glove style. As it is impossible to test every situation encountered and all chemical and mixtures of chemicals, the user is advised to test the particular glove style before and periodically during use.